



**ZYMO RESEARCH**

*The Beauty of Science is to Make Things Simple*

# INSTRUCTION MANUAL

## **Quick-gDNA™ Blood MiniPrep**

Catalog Nos. **D3072 & D3073**

### **Highlights**

- Quick purification of high quality DNA from *whole blood, plasma, and serum* in less than 15 minutes using innovative *Fast-Spin* column technology.
- Compatible with commonly used anticoagulants (i.e., EDTA, heparin, citrate).
- Unique extraction technology excludes the use of Proteinase K and organic denaturants.
- Isolated DNA is ideal for PCR, endonuclease digestion, bisulfite conversion/methylation detection, sequencing, genotyping, etc.

### **Contents**

Product Contents .....	1
Specifications.....	1
Product Description.....	2
Buffer Preparation.....	3
Protocol.....	3
Troubleshooting .....	4
Ordering Information .....	4
Related Products .....	5

Satisfaction of all Zymo Research products is guaranteed. If you should be dissatisfied with this product please call 1-888-882-9682.

Use the **Quick-gDNA™ Blood MicroPrep** (D3070, D3071) for the recovery of up to 5 µg/prep DNA from small samples.

For high-throughput purification (96-well, 5 µg DNA/well) use the **ZR-96 Quick-gDNA™ Blood** (D3075, D3076, D3077).

## Product Contents

<b>Quick-gDNA™ Blood MiniPrep (Kit Size)</b>	<b>D3072 (50 preps.)</b>	<b>D3073 (200 preps.)</b>	<b>Storage Temperature</b>
<b>Genomic Lysis Buffer*</b>	50 ml	2x 100ml	Room Temp.
<b>DNA Pre-Wash Buffer**</b>	15 ml	50 ml	Room Temp.
<b>g-DNA Wash Buffer</b>	50 ml	100 ml	Room Temp.
<b>DNA Elution Buffer</b>	10 ml	2x 10 ml	Room Temp.
<b>Zymo-Spin™ IIC Columns</b>	50	200	Room Temp.
<b>Collection Tubes</b>	100	400	Room Temp.
<b>Instruction Manual</b>	1	1	-

Note - Integrity of kit components is guaranteed for up to one year from date of purchase. Reagents are routinely tested on a lot-to-lot basis to ensure they provide the highest performance and reliability.

\* *Recommended:* Add beta-mercaptoethanol to 0.5%(v/v) i.e., 250 µl per 50 ml or 500 µl per 100 ml.

\*\* A precipitate may have formed in the DNA Pre-Wash Buffer during shipping. To completely resuspend the buffer, incubate the bottle at 30-37 °C for 30 minutes and mix by inversion. DO NOT MICROWAVE.

## Specifications

- **Sample Sources** – Whole blood, plasma, or serum from humans, mice, rats, etc.
- **DNA Purity** – High-quality DNA is eluted with **DNA Elution Buffer** or water. DNA is especially well suited for PCR and other downstream applications.  $A_{260}/A_{280} > 1.8$
- **Size Limits** – Capable of recovering genomic DNA up to and above 40 kb. In most instances, mitochondrial DNA and viral DNA (if present) will also be recovered.
- **DNA Recovery** – Up to 25 µg total DNA is eluted into ≥50 µl (30 µl minimum) **DNA Elution Buffer** or water. Human whole blood will typically yield 3-7 µg DNA per 100 µl blood sampled.
- **Product Detergent Tolerance** – ≤5% Triton X-100, ≤5% Tween-20, ≤5% Sarkosyl, ≤0.1% SDS.
- **Equipment Needed** – Microcentrifuge

Note - ™ Trademarks of Zymo Research Corporation. This product is for research use only and should only be used by trained professionals. It is not intended for use in diagnostic procedures. Some reagents included with this kit are irritants. Wear protective gloves and eye protection. Follow the safety guidelines and rules enacted by your research institution or facility.

### ZYMO RESEARCH CORP.

Phone: (949) 679-1190 ▪ Toll Free: (888) 882-9682 ▪ Fax: (949) 266-9452 ▪ info@zymoresearch.com ▪ www.zymoresearch.com

## Product Description


The **Quick-gDNA™ Blood MiniPrep** is a simple procedure for the rapid isolation of total DNA (e.g., genomic, mitochondrial, viral) from a variety of biological sample sources. This product has been optimized for maximal recovery of ultra-pure DNA without RNA contamination and is compatible with whole blood (fresh or stored), serum, and plasma.

For processing, simply add the specially formulated **Genomic Lysis Buffer** to a sample, vortex, and transfer the mixture to the supplied **Zymo-Spin™ Column**. There is no need for organic denaturants or Proteinase K digestion because of the unique chemistries featured in the kit. Instead, the product features *Fast-Spin* technology to yield high-quality, purified DNA in just minutes (see below). PCR inhibitors are effectively removed during the purification process. DNA purified using the **Quick-gDNA™ Blood MiniPrep** is suitable for PCR, nucleotide blotting, DNA sequencing, restriction endonuclease digestion, bisulfite conversion/methylation analysis, and other downstream applications.



### Ultra-pure DNA for...

- ✓ PCR
- ✓ Endonuclease Digestion
- ✓ Genotyping
- ✓ Bisulfite Conversion  
& Methylation Analysis

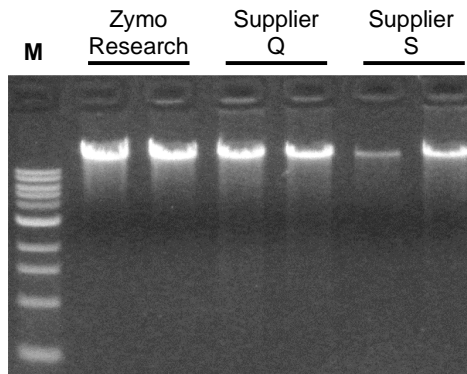

High yield/quality DNA is successfully isolated from porcine whole blood using the **Quick-gDNA™ Blood MiniPrep** (D3072). Equivalent amounts (100 µl) of blood were processed without Proteinase K using the **Quick-gDNA™ Blood MiniPrep** in half the time as compared to the kits from suppliers Q and S. Equal volumes of eluted DNA were then analyzed (in duplicate) in a 0.8% (w/v) TAE/agarose/ethidium bromide gel. The size marker "M" is a 1 kb ladder (Zymo

The **ZymoBead™ Genomic DNA Kit (D3004, D3005)** and **ZR Serum DNA Kit™ (D3013)** are recommended for scaleable DNA yields >25 µg/prep. Both feature silica beads instead of a spin column.

Zymo Research offers the following for rapid, precise DNA methylation detection...

- 1.) **EZ DNA Methylation™ Kit** (D5001, D5002, D5003, D5004)
- 2.) **EZ DNA Methylation-Gold™ Kit** (D5005, D5006, D5007, D5008)
- 3.) **EZ DNA Methylation-Direct™ Kit** (D5020, D5021, D5022, D5023)

## ZYMO RESEARCH CORP.

Phone: (949) 679-1190 ▪ Toll Free: (888) 882-9682 ▪ Fax: (949) 266-9452 ▪ info@zymoresearch.com ▪ www.zymoresearch.com

For **Technical Assistance**,  
please contact 1-888-882-9682  
or E-mail  
tech@zymoresearch.com.

## **Buffer Preparation**

- ✓ **Recommended:** Add beta-mercaptoethanol (user supplied) to the **Genomic Lysis Buffer** to a final dilution of 0.5%(v/v) i.e., 250 µl per 50 ml or 500 µl per 100 ml.

## **Protocol**

*The following is for the purification of DNA from 100 µl whole blood, serum or plasma (the volumes can be adjusted up to 200 µl (max.) depending on your requirements). Fresh, frozen, or preserved blood (in EDTA, citrate, or heparin) can be used. If material cannot be processed immediately, the sample can be “stabilized” for later processing (as noted below) although the immediate processing of blood samples is recommended.*

1. Add 400 µl of **Genomic Lysis Buffer** to 100 µl of blood, serum, or plasma (4:1). Mix completely by vortexing 4-6 seconds, then let stand 5-10 minutes at room temperature.<sup>1</sup>

**Note:** Add 200 µl Genomic Lysis Buffer to all samples <50 µl. For samples larger than 50 µl, add a proportional amount (4:1) of Genomic Lysis Buffer (e.g., Add 800 µl Genomic Lysis Buffer to 200 µl blood).

2. Transfer the mixture to a **Zymo-Spin IIC™ Column<sup>2</sup>** in a **Collection Tube**. Centrifuge at 10,000 x g for one minute. Discard the **Collection Tube** with the flow through.
3. Transfer the **Zymo-Spin™ IIC Column** to a new **Collection Tube**. Add 200 µl of **DNA Pre-Wash Buffer** to the spin column. Centrifuge at 10,000 x g for one minute.
4. Add 500 µl of **g-DNA Wash Buffer** to the spin column. Centrifuge at 10,000 x g for one minute.
5. Transfer the spin column to a clean microcentrifuge tube. Add ≥50 µl **DNA Elution Buffer** or water to the spin column<sup>3</sup>. Incubate 2-5 minutes at room temperature and then centrifuge at top speed for 30 seconds to elute the DNA. The eluted DNA can be used immediately for molecular based applications or stored ≤-20°C for future use.

### **Notes:**

<sup>1</sup> For the inclusion of **small DNAs** from serum, add 0.3 volumes isopropanol to the mixture. (For example, to a 800 µl mixture of serum and Genomic Lysis Buffer add 240 µl isopropanol.)

<sup>2</sup> The column capacity is ~1 ml.

<sup>3</sup> Elution of DNA from the column is dependent on pH and temperature. If water is used, ensure the pH is >6.0. Also, the total yield may be improved by eluting the DNA with Elution Buffer or water pre-equilibrated to 60-70°C.

**Delayed Processing (Stabilization) of Blood Samples:** The immediate processing of blood with this kit is recommended. However, if blood cannot be processed immediately, samples can be “stabilized” in **Genomic Lysis Buffer** for processing at a later time. To do this, add *four* volumes of **Genomic Lysis Buffer** to *each* volume of whole blood (4:1), then vortex. Blood samples mixed with **Genomic Lysis Buffer** can be stored at room temperature for 1-2 weeks, 0-4°C for 1-2 months, -20°C for 6 months to a year, or <-70°C for many years. Samples stored at ≤4°C should reach room temperature prior to processing. Begin at Step 2 in the standard protocol (above) when purifying DNA from blood samples stabilized in **Genomic Lysis Buffer**.

### **ZYMO RESEARCH CORP.**

Phone: (949) 679-1190 ▪ Toll Free: (888) 882-9682 ▪ Fax: (949) 266-9452 ▪ info@zymoresearch.com ▪ www.zymoresearch.com

## Troubleshooting:

1. **DNA degradation:** Check for DNase contamination. All reagents supplied with the **Quick-gDNA™ Blood MiniPrep** are DNase-free. However, DNase contamination could result during the processing of some samples. Check pipets, pipet tips, microcentrifuge tubes, etc., and exercise the appropriate precautions during the DNA purification procedure.
2. **DNA is not performing well in subsequent experiments:** Ensure the correct volume of **Genomic Lysis Buffer** has been added to the sample. Also, make sure all centrifugation steps are completed for the indicated times and speeds (rcfs). Failure to do so may result in incomplete washing, which may cause salts to be eluted with the DNA affecting quantitation and subsequent experiments including enzymatic processes like PCR.
3. **RNA contamination:** The buffers in this kit are designed to efficiently hydrolyze and remove RNA during the DNA purification procedure.
4. **Avian Blood:** Reduce input to 10-25 µl when working with avian blood. Add 7 volumes of **Genomic Lysis Buffer** to the sample. Thoroughly mix the sample by vortexing intermittently for 5 minutes. Transfer to the **Zymo-Spin™** column and proceed with step 2 of the protocol.

## Ordering Information

Product Description	Cat. No.	Kit Size
<b>Quick-gDNA™ Blood MicroPrep</b>	D3070	50 preps.
	D3071	200 preps.
<b>Quick-gDNA™ Blood MiniPrep</b>	D3072	50 preps.
	D3073	200 preps.
<b>Quick-gDNA™ Blood MidiPrep</b>	D3074	25 preps.
<b>ZR-96 Quick-gDNA™ Blood</b>	D3075	2x 96 well
	D3076	4x 96 well
	D3077	10x 96 well

For Individual Sale	Cat. No.	Amount
<b>Genomic Lysis Buffer</b>	D3004-1-50	50 ml
	D3004-1-100	100 ml
<b>DNA Pre-Wash Buffer</b>	D3004-5-15	15 ml
	D3004-5-30	30 ml
	D3004-5-50	50 ml
<b>g-DNA Wash Buffer</b>	D3004-2-50	50 ml
	D3004-2-100	100 ml
<b>DNA Elution Buffer</b>	D3004-4-10	10 ml
<b>Zymo-Spin™ IIC Columns</b>	C1011-50	50
	C1011-250	250
<b>Collection Tubes</b>	C1001-50	50
	C1001-500	500
	C1001-1000	1,000

### ZYMO RESEARCH CORP.

## Popular Products From Zymo Research

Product	Description	Kit Size (Preps)	Catalog No. (Format)
<b>Fragment DNA Purification</b>			
<b>DNA Clean &amp; Concentrator™-5</b>	Clean and concentrate up to 5µg DNA into ≥6 µl elution volume in as little as 2 minutes with no wash residue carryover.	50 200 50 200	<b>D4003</b> (uncapped) <b>D4004</b> (uncapped) <b>D4013</b> (capped) <b>D4014</b> (capped)
<b>DNA Clean &amp; Concentrator™-25</b>	Clean & concentrate 25 µg of DNA into ≥25 µl elution volume in as little as 2 minutes with no wash residue carryover.	50 200 50 200	<b>D4005</b> (uncapped) <b>D4006</b> (uncapped) <b>D4033</b> (capped) <b>D4034</b> (capped)
<b>ZR-96 DNA Clean &amp; Concentrator™-5</b>	Quick (15 minute), high-output recovery of up to 5 µg pure DNA into 10-15 µl minimum elution volume allows for highly concentrated DNA.	2x96 4x96	<b>D4023</b> <b>D4024</b>
<b>Genomic DNA Clean &amp; Concentrator™</b>	Quick (5 minute) clean-up of up to 10 µg high molecular weight DNA (≤200 kb) from any enzymatic reaction or impure preparation without precipitations.	25 100	<b>D4010</b> (capped) <b>D4011</b> (capped)
<b>Zymoclean™ Gel DNA Recovery Kit</b>	Purify DNA from high and low-melting agarose gels in minutes	50 200 50 200	<b>D4001</b> (uncapped) <b>D4002</b> (uncapped) <b>D4007</b> (capped) <b>D4008</b> (capped)
<b>ZR-96 Zymoclean™ Gel DNA Recovery Kit</b>	High-throughput DNA purification from high and low-melting agarose gels.	2x96 4x96	<b>D4021</b> <b>D4022</b>
<b>Zymoclean™ Large Fragment DNA Recovery Kit</b>	Purify high molecular weight DNA (≤200 kb) from high and low-melting agarose gels in minutes	25 100	<b>D4045</b> (capped) <b>D4046</b> (capped)
<b>OneStep™ PCR Inhibitor Removal Kit</b>	Fast, one step procedure for removal of PCR inhibitors such as polyphenolics, humic/fulvic acids, melanin, etc. for successful PCR and other downstream applications.	50 2x96	<b>D6030</b> <b>D6035</b>
<b>Plasmid DNA Purification</b>			
<b>Zyppy™ Plasmid Miniprep Kit</b>	Pellet-Free™ plasmid DNA purification in less than 10 minutes. Recover up to 25 µg DNA in as low as 30 µl.	50 100 400	<b>D4036</b> <b>D4019</b> <b>D4020</b>
<b>Zyppy™-96 Plasmid Miniprep</b>	The fastest and simplest high-throughput method for plasmid purification.	2x96 4x96 8x96	<b>D4041</b> <b>D4042</b> <b>D4043</b>
<b>Zyppy™ Plasmid Midiprep Kit</b>	Pellet-Free™ plasmid DNA purification in 15 minutes in a 150 µl minimum elution volume).	25 50	<b>D4025</b> <b>D4026</b>
<b>ZR Plasmid MiniPrep™ Classic</b>	Plasmid DNA purification in minutes: (alkaline lysis/spin column format for low 30 µl elution volume).	50 100 400	<b>D4036</b> <b>D4019</b> <b>D4020</b>
<b>Genomic DNA Purification</b>			
<b>Quick-gDNA™ MiniPrep</b>	Easy purification of genomic DNA from whole blood, plasma, serum, body fluids, buffy coat, lymphocytes, tissue, swabs or cultured cells in as little as 15 minutes <u>without</u> the use of Proteinase K or organic denaturants.	50 200 50 200	<b>D3006</b> (uncapped) <b>D3007</b> (uncapped) <b>D3024</b> (capped) <b>D3025</b> (capped)
<b>ZR-96 Quick-gDNA™</b>	Simple, high throughput (96-well) purification of DNA from whole blood, plasma, serum, body fluids, buffy coat, lymphocytes, tissue, swabs, or cultured cells in about 30 minutes.	2x96 4x96 10x96.	<b>D3010</b> <b>D3011</b> <b>D3012</b>
<b>ZR-Genomic DNA™-Tissue MiniPrep</b>	For high quality DNA purification from <u>solid tissues</u> (e.g., tail snips, ear punches, adipose tissue, etc.), body fluids, cultured cells, buccal cells, FFPE tissues, hair, and other biological sources using Proteinase K and Fast.	50 200	<b>D3050</b> <b>D3051</b>
<b>Environmental DNA Purification Kits</b>	Unique BashingBead™ technology allows isolation of DNA from samples refractory to conventional lysis procedures including tough-to-lyse tissues, soil samples, feces, plants, seeds, insects, bacteria, yeast, filamentous fungi, unicellular and filamentous algae, and protozoa		<b>Visit website for a comprehensive list</b>

Please visit our website to see our complete line-up of products.

### ZYMO RESEARCH CORP.

Phone: (949) 679-1190 ▪ Toll Free: (888) 882-9682 ▪ Fax: (949) 266-9452 ▪ info@zymoresearch.com ▪ www.zymoresearch.com